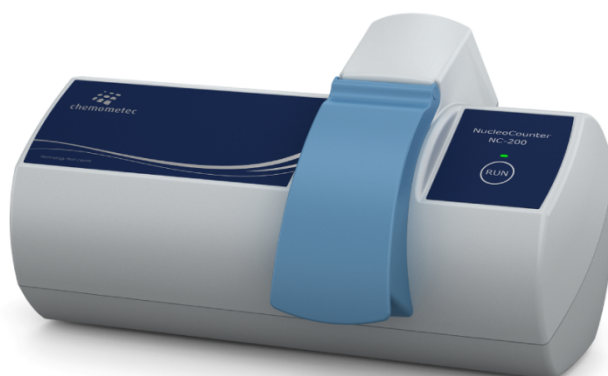


# NucleoCounter® Lids

Package insert for 929-0007, 929-0008, 929-0009 and 929-0010.



**Figure 1. NucleoCounter with large lid for storage.**

The NucleoCounter instrument is delivered with two lids. The larger of the two lids, shown on the figure above, protects the insertion area from dust and dirt. This lid should be placed on the instrument in the closed position when the instrument is not in use. The smaller of the two lids (see the figure below) protects the insertion area against ambient light. This lid is easier to handle and can replace the large lid if it is more convenient during a measurement session. However, for protection replace the small lid with the large lid, when the session is over.



**Figure 2. NucleoCounter with the small version of the lid, which can be used when measuring samples.**